


# Download Instructions for Thermo Scientific SkanIt Software Protocols of the Thermo Scientific Multiskan GO

## 1) Downloading the protocols from the website

1. Click the protocol download icon and choose **Save**.

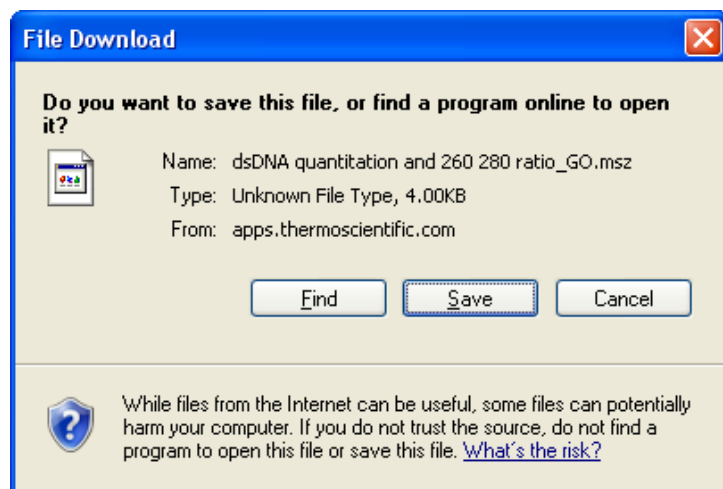

2. Save the file on your computer.

## 2) Importing the downloaded protocols to the SkanIt Software database

1. Open SkanIt Software for Multiskan GO.
2. Select **Import** from the **Home** view:



3. Search the protocol file by clicking **Browse**. Select the protocol file (\*\_GO.msz) and click **Open**.
4. Click **Next** to continue.
5. Select the SkanIt Software database folder where you want the protocol to be imported to. Click **Finish**. The software will notify when the protocol has been imported successfully.

6. The protocol can now be opened either by clicking **Open** and searching for the imported protocol, or by clicking the protocol name from the **Recent Sessions** list in the **Home** view.

7. Before starting the measurement:

- Check that the pipetting order of the samples on the microplate matches the sample definitions in the **Layout** view.
- Check that the measurement wavelength in the **Protocol** view is correct.
- Check that the calculations in the **Results** view are correct.